



La Pias Ear Tag Application

- Using the applicator, detach the female portion of the ear tag and place it into the red slot of the applicator.
 - It is recommended to clean the tags and applicator with 70% alcohol or UV irradiation before application to prevent infection.



- After making sure the female portion is properly inserted, place the male portion of the tag into the appropriate slot of the applicator (opposite the female).
 - Using forceps may help with placement.



- Avoiding any blood vessels or the outer ring of ear cartilage, place the tag in the center of the ear and apply.
 - Check to be sure the point of the tag has come through the ear and is properly seated in the backing.

