


Waste Gas Exposure Results

Personal dosimeter badges were worn during anesthesia via face mask at 2.5% isoflurane for 30 minutes. Vaporizers were operated at manufacturer recommended flow rate settings — SomnoSuite: 27±2 mL/min (1.5x animal's minute volume as determined by animal body weight); Traditional Vaporizer: 500 mL/min.

- SomnoSuite resulted in undetectable exposure — below minimum dosimeter limits (n=3)
- Traditional vaporizer resulted in exposure levels above NIOSH guidelines (n=3)

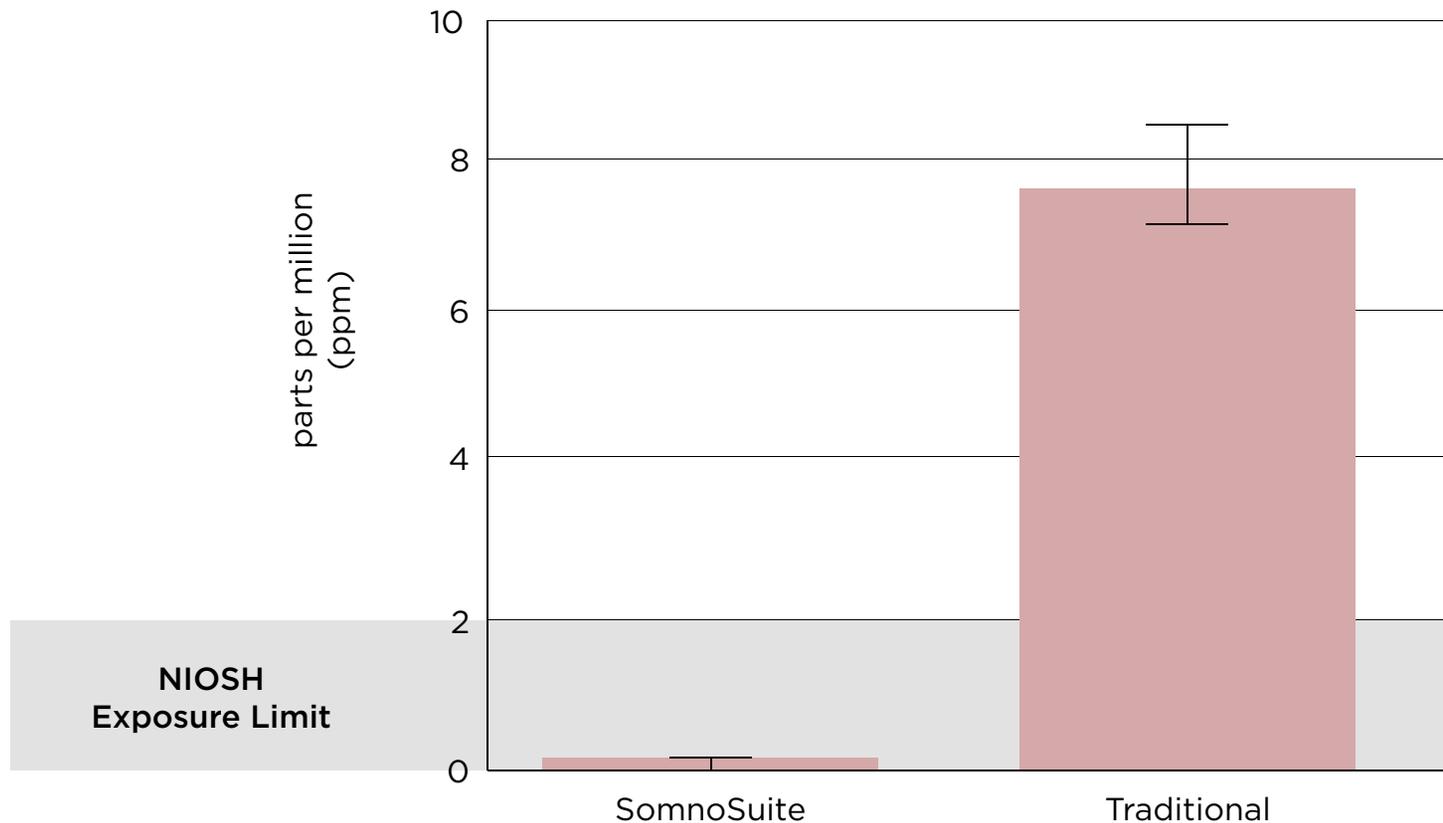

For more information contact us:

888.572.8887

www.kentscientific.com

